


2,4-D CONCISE FLOWCHART

1.



Add 500 μ L of either Standards (calibrators), Control or Samples to the bottom of each test tube.

5.

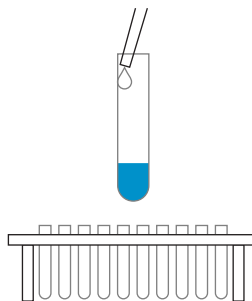

Add 500 μ L of Color Reagent down the inside wall of each test tube.

2.



Add 200 μ L of 2,4-D Enzyme Conjugate to each test tube.

Vortex or swirl.

6.



Incubate for 20 minutes.


Prepare blank.

3.



Incubate for 20 minutes.

7.

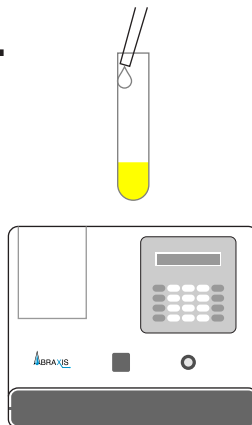

Add 500 μ L of Stopping Solution to each test tube.

Read OD 450

4.



Flood tubes with deionized water.

Invert tubes and blot.

Repeat this step 4 times.

For Ordering or Technical Assistance Contact:
Eurofins Abraxis
124 Railroad Drive, Warminster, PA 18974 USA
Phone: 215-357-3911 Fax: 215-357-5232
E-mail: info.ET.Warminster@eurofinsUS.com
www.abraxiskits.com

2,4-D Tube Kit Part # 54004B, 40 Test